


MESTRA®
CALENTADOR DE PLACAS BASE
WATER BATH
RÉCHAUFFEUR DE PLAQUES BASE
R-080078

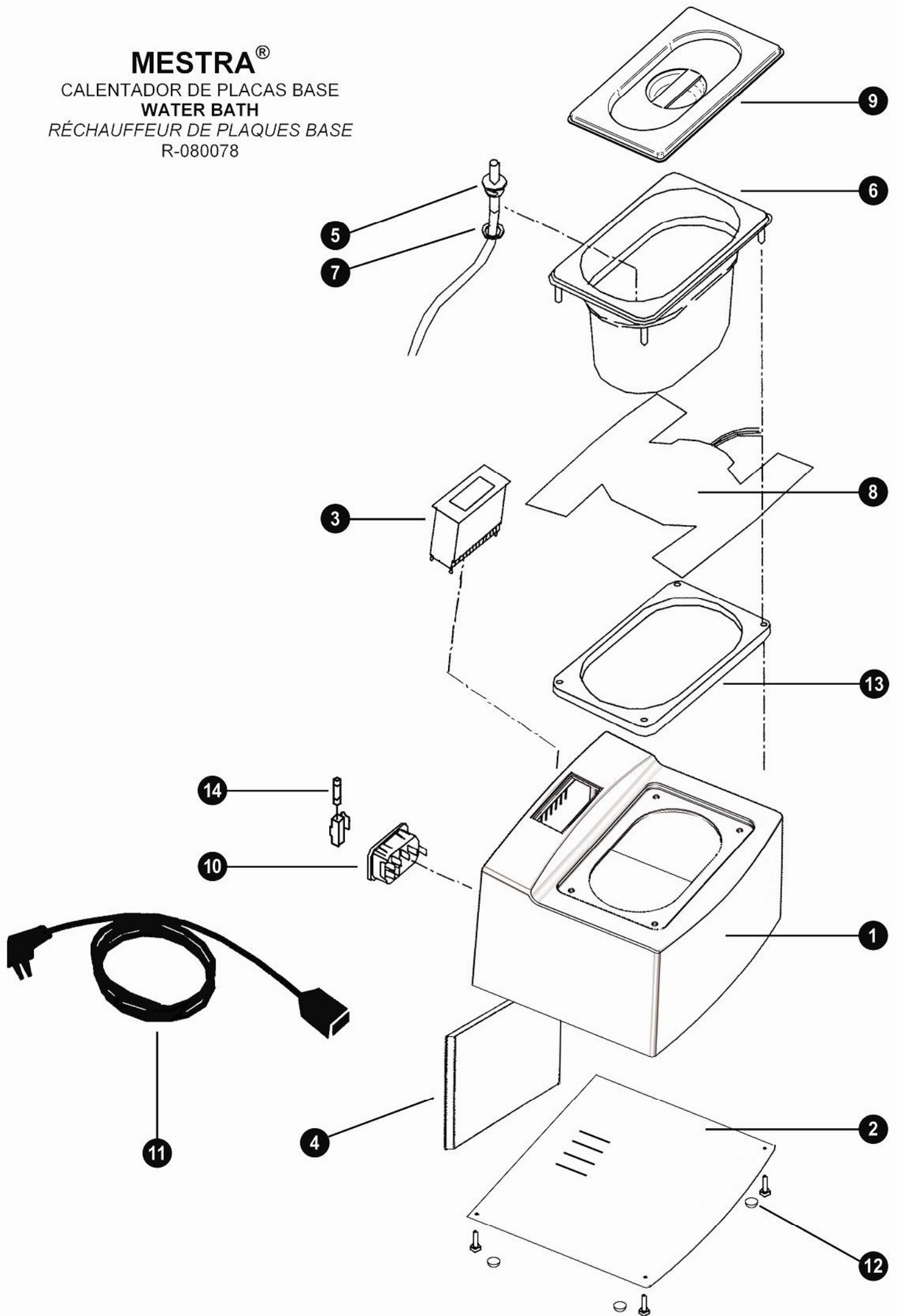

MESTRA®
 CALENTADOR DE PLACAS BASE
WATER BATH
 RÉCHAUFFEUR DE PLAQUES BASE
 R-080078

Nº	Ref.	Descripción	Description
1	080078-01	Carcasa	Housing
2	080078-02	Tapa inferior	Lower lid
3	080078-03	Control 230V	Control 230V
4	080078-04	Fibra aislante	Insulation fiber
5	080078-05	Sonda	Termal probe
6	080077-01	Cuba	Container
7	040210-07	Junta tórica	O-ring
8	080077-04	Resistencia 230V	Resistor 230V
9	080077-09	Tapa superior	Upper lid
10	080520-16	Base de enchufe	Connection base
11	000250	Cable 230V	Cable 230V
12	050364-04	Pata	Leg
13	080077-05	Junta de silicona	Silicone gasket
14	000231	Fusible 4A	Fuse 4A